


MycoFog™ Quick Start guide

Please consult full manual at <https://mycofog.com>

Day before use: Charge MycoFog™ Device with supplied cable.

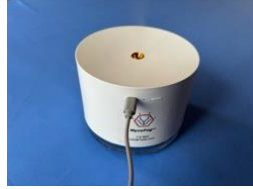


> 8 hours

Day of use: Take out all cell cultures from incubator, set Incubator to **37°C** or **45°C**, turn off gas control and remove water pan or turn off active humidification.

1. Open MycoFog™



2. Pour MycoFog™ Reagent into reagent reservoir of MycoFog™

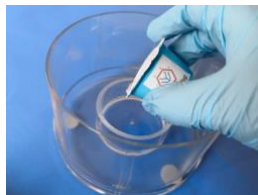

3. Prime wick for 60 seconds in reagent reservoir of MycoFog™



4. Insert primed end of wick into upper compartment by removing and replacing protector over wick, press firmly. Replace top of MycoFog™.



5. Place MycoFog™ in incubator, mid-shelf, wait 5 minutes.



5 minutes

6. Press Start



7. Close incubator door & set timer.



180 Minutes (or **240** mins if using 37 °C)

8. Remove MycoFog™, *if necessary* return temperature to 37°C. Your Incubator is now decontaminated and ready for use!